

## PEEK & Polymer Chemical Resistance

### PEEK & Polymer Chemical Resistance

PEEK exhibits excellent resistance to a wide range of organic and inorganic chemicals. The compatibility of PEEK with many chemicals at 20°C (68°F) has been investigated and the results for unreinforced grades are given in the table below.

PEEK is compatible with almost any of the solvents used in HPLC. The only solvent which will attack PEEK are

concentrated nitric acid and sulphuric acids. However, PEEK tubing can safely withstand 20–30% nitric acid when passivating a system.

Methylene Chloride, DMSO and THF may cause swelling in PEEK. The highest temperature we recommend for PEEK is 100°C. Up to this temperature the tubing will maintain the pressure rating stated.

A = suitable

B = marginal-dependent on application

C = not recommended

#### PEEK and Polymer chemical resistance

Resistance at 20°C	PEEK	PE	PP	PPS	FEP	ETFE	PFA	PTFE
Acetaldehyde	A				A		A	
Acetic acid (20%)	A	A	A	A	A	A	A	A
Acetic acid (80%)	A	A	A	A	A		A	A
Acetic acid (glacial)	A	A	A	A	A	A	A	A
Acetone	A	B	A	A	A	A	A	A
Acetonitrile	A			A	A	A	A	A
Acrylic acid	A			A	A		A	
Ammonia, anhydrous	A			A	A		A	
Ammonia (10%)	A	B	A	A	A	A	A	A
Ammonia (Liquid)	B			A	A		A	
Ammonium hydroxide	A	A	A	A	A	A	A	A
Aqua regia	C			C	A		A	
Aromatic hydrocarbons	A	B	C	A	A		A	
Benzene	A	B	C	A	A	A	A	A
Benzoic acid	A			A	A		A	
Benzaldehyde	A			A	A		A	
Bromine/dibromoethane	C			C	A		A	
Bromine (dry)	C			C	A		A	
Bromine (wet)	C			C	A		A	
Boric acid	A			A	A		A	
Butanol	A	A	A	A	A	A	A	A
Calcium hydroxide	A			A	A		A	
Carbon tetrachloride	A			A	A		A	
Chlorine (gas)	A			C	A		A	
Chlorine (liquid)	C			C	A		A	
Chloroacetic acid	A	B	B	A	A	A	A	A
Chlorobenzene	A			A	A		A	
Chloroform	A	B	B	A	A	A	A	A
Cyclohexane	A	B	C	A	A	A	A	A
Cyclohexanone	A	C	C	A	A	A	A	A
Diethylamine	A	C	A		A	A	A	A
Diethylether	A			A	A		A	
Diethylformamide	A	A	A	A	A	A	A	A
Dioxane	A			A	A	A	A	
Ethanol	A	B	A	A	A	A	A	A
Ether	A	B	C	A	A	A	A	A
Ethyl acetate	A	B	A	A	A	A	A	A
Ethylene chloride		B	B	A	A	A	A	A
Ethylene glycol	A	A	A	A	A	A	A	A
Heptane	A	B	B	A	A	A	A	A

## PEEK &amp; Polymer Chemical Resistance

Resistance at 20 °C	PEEK	PE	PP	PPS	FEP	ETFE	PFA	PTFE
Hexane	A	B	B	A	A	A	A	A
Hydrobromic acid (100%)	C	B	B	A	A		A	A
Hydrobromic acid (20%)	C	B	A		A	A	A	A
Hydrochloric acid (100%)	A		B	C	A	A	A	A
Hydrochloric acid (20%)	A	A	B	C	A	A	A	A
Hydrofluoric acid (100%)	C		B	C	A	A	A	A
Hydrofluoric acid (20%)		A	A	A	A	A	A	A
Hydrogen peroxide (100%)	A	B	B	C	A		A	A
Hydrogen peroxide (50%)	A	B	B		A	A	A	A
Hydrogen peroxide (10%)	A	A	A	A	A	A	A	A
Iso-octane	A		A	A	A		A	A
Isopropanol	A	A	A	A	A		A	A
Isopropyl ether		A	B		A		A	A
Ketones	A	B	B	A	A		A	A
Methanol	A	A	A	A	A	A	A	A
Methyl dichloride			C		A		A	
Methyl ethylketone	A	B	B	A	A	A	A	A
Methylene chloride	B	B	B	A	A	A	A	A
Nitric acid (100%)	C	B	C	C	A	A	A	A
Nitric acid (20%)	A	B	A	C	A	A	A	A
Pentane	A	C	C		A		A	A
Perchloric acid	A	B	B		B		B	A
Phenol (dilute)	A			A	A		A	
Phenol (conc.)	C			A	A		A	
Phosphoric acid (100%)	A	B	A	A	A	A	A	A
Phosphoric acid (40%)	A	A	A	A	A	A	A	A
Potassium hydroxide (dilute)	A			A	A		A	
Potassium hydroxide (70%)	A			A	A		A	
Propanol	A				A		A	
Pyridine	A	B	A	A	A	A	A	A
Sodium hydroxide (80%)	A	B	A	A	A		A	A
Sodium hydroxide (20%)	A	A	A	A	A	A	A	A
Sulphuric acid (100%)	C	B	B	A	A	A	A	A
Sulphuric acid (75%)	C	A	A	A	A	A	A	A
Sulphuric acid (40%)	A	A	A	A	A	A	A	A
Tetrahydrofuran	B	B	B	A	A	A	A	A
Toluene	A	B	B	A	A	A	A	A
Trichloroacetic acid	A	A	A	A	A	A	A	A
Trichloroethane	A		B		A		A	A
Trichloroethylene	A			A	A		A	
Triethylamine			C		A	A	A	A
Water	A	A	A	A	A	A	A	A
Water (distilled)	A	A	A	A	A	A	A	A
Xylene	A			A	A		A	

## The chemical resistance of PEEK at high temperature

Chemical	Temperature	Rating	Chemical	Temperature	Rating
Acetic acid	200 °C (420 °F)	B	Methane	200 °C (420 °F)	A
Ammonia gas	200 °C (420 °F)	A	Methylethylketone	200 °C (420 °F)	C
Carbon monoxide gas	200 °C (420 °F)	A	Nitrobenzene	200 °C (420 °F)	C
Ethylene glycol	200 °C (420 °F)	B	Phosphoric acid (50%)	200 °C (420 °F)	A
Ethylene glycol (50%)	140 °C (284 °F)	A	Sodium hydroxide solution	200 °C (420 °F)	A
Hydrogen sulphide gas	200 °C (420 °F)	A	Sulphuric acid (50%)	200 °C (420 °F)	B
Liquid ammonia	200 °C (420 °F)	B	Sulphur dioxide gas	200 °C (420 °F)	A